

## Human MCP-1 ELISA Kit

## EH2MCP1 EH2MCP12

1572.3

<b>Number</b>	<b>Description</b>
<b>EH2MCP1</b>	<b>Human MCP-1 ELISA Kit</b> , sufficient reagents for 96 determinations <b>Kit Contents:</b> <b>Anti-Human MCP-1 Precoated 96-well Strip Plate</b> , 1 each <b>Lyophilized Recombinant Human MCP-1 Standard</b> , 2 vials <b>Standard Diluent</b> , 25 ml, contains 0.1% sodium azide <b>Biotinylated Antibody Reagent</b> , 12 ml, contains 0.1% sodium azide <b>30X Wash Buffer</b> , 50 ml <b>Streptavidin-HRP Concentrate</b> , 75 µl <b>Streptavidin-HRP Dilution Buffer</b> , 14 ml <b>TMB Substrate</b> , 13 ml <b>Stop Solution</b> , 13 ml, contains 0.16 M sulfuric acid <b>Adhesive Plate Covers</b> , 6 each
<b>EH2MCP12</b>	<b>Human MCP-1 ELISA Kit</b> , sufficient reagents for 2 × 96 determinations <b>Kit Contents:</b> <b>Anti-Human MCP-1 Precoated 96-well Strip Plate</b> , 2 each <b>Lyophilized Recombinant Human MCP-1 Standard</b> , 4 vials <b>Standard Diluent</b> , 2 × 25 ml, contains 0.1% sodium azide <b>Biotinylated Antibody Reagent</b> , 2 × 12 ml, contains 0.1% sodium azide <b>30X Wash Buffer</b> , 2 × 50 ml <b>Streptavidin-HRP Concentrate</b> , 2 × 75 µl <b>Streptavidin-HRP Dilution Buffer</b> , 2 × 14 ml <b>TMB Substrate</b> , 2 × 13 ml <b>Stop Solution</b> , 2 × 13 ml, contains 0.16 M sulfuric acid <b>Adhesive Plate Covers</b> , 12 each

For research use only – not for use in diagnostic procedures.

**Storage:** For maximum stability, store in a non-defrosting -20°C freezer and refer to the expiration date for frozen storage on the label. Alternatively, store at 2-8°C and refer to the expiration date for refrigerated storage. Once thawed, store at 4°C until the expiration date for refrigerated storage. Kit is shipped on dry ice.


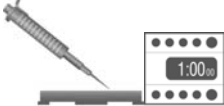

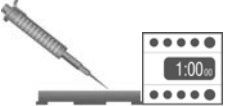
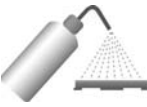

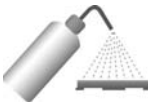





## Introduction

The Thermo Scientific Human Monocyte Chemoattractant Protein-1 (MCP-1) ELISA is an enzyme-linked immunosorbent assay for quantification of human MCP-1 in serum; EDTA and heparin plasma; and culture supernatant.

## Table of Contents

Introduction .....	1
Procedure Summary.....	2
Additional Materials Required.....	2
Precautions.....	3
Sample Preparation.....	3
Reagent Preparation.....	4
Assay Procedure .....	5
Performance Characteristics .....	6
Reference .....	7
Data Templates .....	8

## Procedure Summary

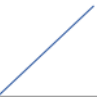- |  |  |   |   |
|--|--|---|---|
| <br><b>1.</b> Add 50 µl of Standard Diluent to each well.             | <br><b>2.</b> Add 50 µl of standards or samples in duplicate. Cover plate and incubate at room temperature (20-25°C) for 1 hour.          | <br><b>3.</b> Wash the plate THREE times.                                       | <br><b>4.</b> Add 100 µl of Biotinylated Antibody Reagent to each well. Cover plate and incubate at room temperature for 1 hour. |
| <br><b>5.</b> Wash the plate THREE times.                             | <br><b>6.</b> Add 100 µl of prepared Streptavidin-HRP Solution to each well. Cover plate and incubate at room temperature for 30 minutes. | <br><b>7.</b> Wash the plate THREE times.                                       | <br><b>8.</b> Add 100 µl Premixed TMB Substrate Solution to each well.   |
| <br><b>9.</b> Develop the plate at room temperature for 20 minutes. | <br><b>10.</b> Add 100 µl of Stop Solution to each well.  | <br><b>11.</b> Measure absorbance. Subtract 550 nm values from 450 nm values. | <br><b>12.</b> Calculate results using curve-fitting statistical software.   |

## Additional Materials Required

- Precision pipettors with disposable plastic tips to deliver 5-1,000 µl and plastic pipettes to deliver 5-15 ml
- A glass or plastic two-liter container to prepare Wash Buffer
- A squirt wash bottle or an automated 96-well plate washer
- 1.5 ml polypropylene or polyethylene tubes to prepare standards – do not use polystyrene, polycarbonate or glass tubes
- Disposable reagent reservoirs
- 15 ml plastic tube to prepare Streptavidin-HRP Solution
- A standard ELISA reader for measuring absorbance at 450 nm and 550 nm. If a 550 nm filter is not available, the absorbance may be measured at 450 nm only. Refer to the instruction manual supplied with the instrument being used.
- Graph paper or a computerized curve-fitting statistical software package

---

## Precautions

- **All samples and reagents must be at room temperature (20-25°C) before use in the assay.**
- Review all instructions carefully and verify components against the Kit Contents list (page 1) before beginning the assay.
- Do not use a heated water bath to thaw samples. Thaw samples at room temperature.
- Sample Diluent must be used when diluting all samples.
- When assaying culture medium, prepare the standard curve and sample dilutions using the same medium used for the cell culture. For example, if RPMI with 10% fetal calf serum (FCS) was used, then use RPMI with 10% FCS to dilute the standard and samples. For best results, use a culture medium that contains a carrier protein such as FCS. The lack of a carrier protein may affect MCP-1 stability and compromise the standard curve. If serum-free media must be used, prepare two standard curves: one with carrier and one without. Compare curves to determine if lack of carrier caused degradation that resulted in a signal decrease.
- If using a multichannel pipettor, always use a new disposable reagent reservoir.
- Use new disposable pipette tips for each transfer and a new adhesive plate cover for each incubation step to avoid cross-contamination.
- Once reagents have been added to the plate, take care NOT to let plate DRY at any time during the assay.
- Avoid microbial contamination of reagents.
- Avoid exposing reagent to excessive heat or light during storage and incubation.
- Do not mix reagents from different kit lots. Discard unused kit components.
- Do not use glass pipettes to measure TMB Substrate Solution. Take care not to contaminate the solution. If the solution is blue before use, DO NOT USE IT.
- Individual components may contain antibiotics and preservatives. Wear gloves while performing the assay to avoid contact with samples and reagents. Please follow proper disposal procedures.
- Some components of this kit contain sodium azide. Please dispose of reagents according to local regulations.

## Sample Preparation

### Sample Handling

- Serum; EDTA or heparin plasma; or culture supernatants may be tested in this assay.
- 50 µl per well of serum, plasma or culture supernatant is required.
- Store samples to be assayed within 24 hours at 2-8°C. For long-term storage, aliquot and freeze samples at -70°C. Avoid repeated freeze-thaw cycles when storing samples.
- Gradually equilibrate samples to room temperature before beginning assay. Do not use a heated water bath to thaw or warm samples. Mix samples by gently inverting tubes.
- If samples are clotted, grossly hemolyzed, lipemic or contaminated, or if there is any question about the integrity of a sample, make a note on the template and interpret results with caution.
- Samples and standards must be assayed in duplicate each time the assay is performed.

### Sample Dilution

- Culture supernatant samples may be evaluated **undiluted** in this assay. If the human MCP-1 concentration possibly exceeds the highest point of the standard curve (i.e., 1,000 pg/ml), prepare one or more 10-fold dilutions of the sample using culture media as the diluent. A 10-fold dilution is prepared by adding 50 µl of sample to 450 µl of culture medium. Mix thoroughly between dilutions.
- Serum and plasma may be evaluated **undiluted** in this assay. If the MCP-1 concentration in a sample possibly exceeds the highest point of the standard curve, (i.e., 1,000 pg/ml), prepare one or more additional five-fold dilutions. A five-fold dilution is prepared by adding 50 µl of sample to 200 µl Standard Diluent. Mix thoroughly between dilutions.

## Reagent Preparation

For procedural differences when using partial plates, look for **(PP)** throughout these instructions.

### Wash Buffer

**Note:** Wash Buffer must be at room temperature before use in the assay. Do not use Wash Buffer if it becomes visibly contaminated during storage.

1. Label a clean glass or plastic two-liter container "Wash Buffer." The 30X Wash Buffer may have a cloudy appearance.
2. If using the 5-plate kit, add 30 ml Wash Buffer to 870 ml water for each plate used, otherwise, add the entire contents of the 30X Wash Buffer bottle (50 ml) to the container and dilute to a final volume of 1.5 L with ultrapure water. Mix thoroughly.

**(PP)** When using partial plates, store the reconstituted Wash Buffer at 2-8°C.


### Standards

- **(PP)** Reconstitute and use one vial of the lyophilized standard per partial plate.
  - Prepare standards just before use and use within one hour of reconstitution. Do not store reconstituted standards.
1. When testing **culture supernatant** samples, reconstitute standard with ultrapure water. Reconstitution volume is stated on the vial label. The standard will dissolve in approximately 1 minute. Mix by gently inverting vial. Use sample culture medium to prepare standard curve serial dilutions.

When testing **serum or plasma** samples, reconstitute standard with ultrapure water. Reconstitution volume is stated on the vial label. Mix by gently inverting the vial until the contents have completely dissolved. Use the Sample Diluent provided to prepare the standard curve serial dilutions.

When testing **serum, plasma or cell culture supernatant samples on the same plate**, validate the medium to establish if the same standard curve can be used for the different sample types. Prepare a standard curve (including a zero/blank) using culture medium to dilute the standard. Use medium containing serum or other protein supplement to maximize stability of the MCP1. Perform this curve in parallel with a standard curve diluted in the Standard Diluent provided. If OD values are within 10% of the mean value for both curves, then the assay may be performed with Standard Diluent, whether testing culture supernatant, plasma, urine or serum samples.

2. Label six tubes, one for each standard curve point: 1,000 pg/ml, 400 pg/ml, 160 pg/ml, 64 pg/ml, 25.6 pg/ml and 0 pg/ml, then prepare serial dilutions for the standard curve as follows:
3. Pipette 330 µl of Standard Diluent into the first tube (i.e., 1,000 pg/ml)
4. Pipette 240 µl of Standard Diluent into the remaining tubes.
5. Pipette 82.5 µl of the reconstituted standard into the first tube (i.e., 1,000 pg/ml) and mix for a 1:5 dilution.
6. Pipette 160 µl of this dilution into the next tube (i.e., 400 pg/ml) and mix for a 1:2.5 dilution.
7. Repeat the 1:2.5 serial dilutions (using 160 µl) three more times to complete the standard curve points.

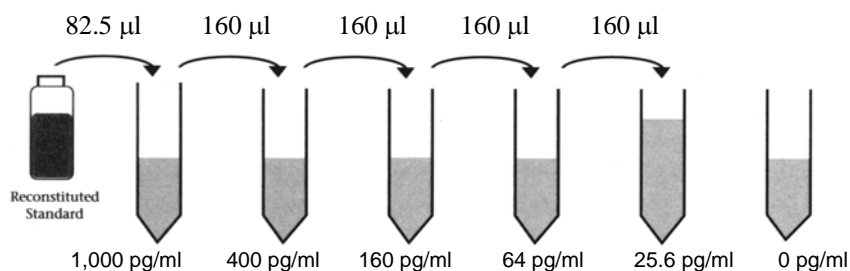

---

## Assay Procedure

### A. Sample Incubation

- **(PP)** Determine the number of strips required and leave these strips in the plate frame. Place remaining unused strips in the provided foil pouch with desiccant and store at 2-8°C. Make sure foil pouch is sealed tightly. After completing assay, retain plate frame for second partial plate. When using the second partial plate, place strips securely in the plate frame.
  - Use the Data Template provided to record locations of the zero standard (blank or negative control), standards and samples. Perform five standard points and one blank in duplicate with each series of unknown samples.
1. Add 50 µl of Standard Diluent to each well.
  2. Add 50 µl of reconstituted standards or samples in duplicate to each well. Mix by gently tapping the plate several times.  
**Note:** If the human MCP-1 concentration in any sample possibly exceeds the highest point on the standard curve, 1,000 pg/ml, see Sample Preparation – Sample Dilution section.
  3. Carefully cover plate with a new adhesive plate cover. Ensure all edges and strips are sealed tightly by running your thumb over edges and down each strip. Incubate for one (1) hour at room temperature, 20-25°C.
  4. Carefully remove adhesive cover, discard plate contents and wash plate as described in the Plate Washing section below.

### B. Plate Washing

1. Gently squeeze the long sides of plate frame before washing to ensure all strips securely remain in the frame.
2. Empty plate contents. Use a squirt wash bottle to vigorously fill each well completely with Wash Buffer, then empty plate contents. Repeat procedure two additional times for a total of THREE washes. Blot plate onto paper towels or other absorbent material.  
**Note:** For automated washing, aspirate all wells and wash THREE times with Wash Buffer, overfilling wells with Wash Buffer. Blot plate onto paper towels or other absorbent material.

### C. Biotinylated Antibody Reagent Incubation

- If using a multichannel pipettor, **use a new reagent reservoir and pipette tips** when adding the Biotinylated Antibody Reagent. Remove from the vial only the amount required for the number of strips being used.
1. Add 100 µl of Biotinylated Antibody Reagent to each well containing sample or standard. Mix well by gently tapping the plate several times.
  2. Carefully attach a new adhesive plate cover. Ensure all edges and strips are tightly sealed by running your thumb over edges and down each strip. Incubate plate for one (1) hour at room temperature, 20-25°C.
  3. Carefully remove the adhesive plate cover and wash THREE times with Wash Buffer as described in the Plate Washing section (section B).

### D. Streptavidin-HRP Solution Preparation and Incubation

- Prepare Streptavidin-HRP Solution **immediately before use**. Do not prepare more Streptavidin-HRP Solution than required. Do not store prepared Streptavidin-HRP Solution.
  - Use a 15 ml plastic tube to prepare Streptavidin-HRP Solution.
  - **Note:** If using a multichannel pipettor, **use new reagent reservoir and pipette tips** when adding the prepared Streptavidin-HRP Solution.
1. Briefly centrifuge the Streptavidin-HRP Concentrate to force entire vial contents to the bottom.
  2. **(PP)** Use only the Streptavidin-HRP Solution amount required for the number of strips being used. For each strip, mix 2.5 µl of Streptavidin-HRP Concentrate with 1 ml of Streptavidin-HRP Dilution Buffer. Store Streptavidin-HRP Concentrate reserved for additional strips at 2-8°C.

For one complete 96-well plate, add 30 µl of Streptavidin-HRP Concentrate to 12 ml of Streptavidin-HRP Dilution Buffer and mix gently.

3. Add 100 µl of prepared Streptavidin-HRP Solution to each well.
4. Carefully attach a new adhesive plate cover, ensuring all edges and strips are tightly sealed. Incubate plate for 30 minutes at room temperature, 20-25°C.
5. Carefully remove the adhesive cover, discard plate contents and wash plate as described in the Plate Washing section.

#### E. Substrate Incubation and Stop Step


- Use new disposable reagent reservoirs when adding TMB Substrate Solution and Stop Solution.
  - Dispense from bottle ONLY amount required, 100 µl per well, for the number of wells being used. Do not use a glass pipette to measure the TMB Substrate Solution.
  - **(PP)** Do not combine leftover substrate with that reserved for the second partial plate. Take care not to contaminate remaining TMB Substrate Solution.
1. Pipette 100 µl of TMB Substrate Solution into each well.
  2. Allow color reaction to develop at room temperature in the dark for 20 minutes. Do not cover plate with aluminum foil or a plate sealer. The substrate reaction yields a blue solution that turns yellow when Stop Solution is added.
  3. After 20 minutes, stop the reaction by adding 100 µl of Stop Solution to each well.

#### F. Absorbance Measurement

1. Evaluate plate within 30 minutes of stopping the reaction. Measure absorbance on an ELISA plate reader set at 450 nm and 550 nm. Subtract 550 nm values from 450 nm values to correct for optical imperfections in the microplate. If 550 nm is not available, measure absorbance at 450 nm only. Omitting the 550 nm measurement results in higher absorbance values.

#### G. Calculation of Results

- Use the standard curve to determine human MCP-1 amount in an unknown sample. Generate the standard curve by plotting the average absorbance obtained for each standard concentration on the vertical (Y) axis vs. the corresponding human MCP-1 concentration (pg/ml) on the horizontal (X) axis.
- Calculate results using graph paper or curve-fitting statistical software with a linear regression or 4-parameter logistic fit. Determine the human MCP-1 amount in each sample by interpolating from the absorbance value (Y axis) to MCP-1 concentration (X axis) using the standard curve.
- If the sample was diluted, multiply the obtained interpolated value by the dilution factor to calculate amount of MCP-1 in the sample.
- Absorbance values obtained for duplicates should be within 10% of the mean value. Carefully consider duplicates with values that differ from the mean by greater than 10%.

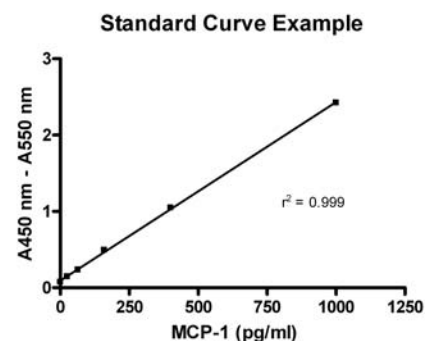

### Performance Characteristics

**Sensitivity:** < 10 pg/ml

The sensitivity or lower limit of detection (LLD)<sup>1</sup> is determined by assaying replicates of zero and the standard curve. The mean signal of zero + 2 standard deviations read in dose from the standard curve is the LLD. This value is the smallest dose that is not zero with 95% confidence.

**Assay Range:** 0-1,000 pg/ml

Suggested standard curve points are 1,000 pg/ml, 400 pg/ml, 160 pg/ml, 64 pg/ml, 25.6 pg/ml, and 0 pg/ml.

#### Reproducibility:

Intra-assay CV: <10%

Inter-assay CV: <10%

**Specificity:** This ELISA is specific for the measurement of natural and recombinant human MCP-1. This ELISA does not cross-react with the following recombinant human chemokines and cytokines: Angiogenin, EGF, FGF-basic, Eotaxin, GM-CSF, GRO $\alpha$ , HB-EGF, HGF, IFN $\alpha$ , IFN $\gamma$ , KGF, MCP-2, MCP-3, MCP-4, MIP-1 $\alpha$ , MIP-1 $\beta$ , PLGF-1, RANTES, TNF $\alpha$ , IL-1 $\alpha$ , IL-1 $\beta$ , IL-1RA, IL-2, IL-3, IL-4, IL-5, IL-6, IL-7, IL-8, IL-9, IL-10, IL-11, IL-12, IL-13, IL-15, IL-16, and IL-17.

**Calibration:** The standards in this ELISA are calibrated to the NIBSC recombinant MCP-1 standard 92/794. One (1) pg of internal standard = 0.58 pg of NIBSC standard = 0.00058 NIBSC units.

**Expected Values:** Serum and plasma samples were collected from apparently healthy human donors and tested in this assay. The levels of human MCP-1 found in each sample type are reported in Table 1.

**Table 1.** Human MCP-1 levels from apparently healthy individuals

<u>Sample Type</u>	<u>Average</u>	<u>Range</u>
Serum (n= 15)	305 pg/ml	228 - 475 pg/ml
EDTA Plasma (n=14)	117 pg/ml	58 - 234 pg/ml
Heparin Plasma (n=14)	278 pg/ml	163 - 438 pg/ml

**Spike and Recovery:** Recovery was determined by spiking various levels of recombinant human MCP-1 into serum and plasma samples collected from apparently healthy individuals and a Standard Diluent control buffer. Mean recoveries are reported in Table 2.

**Table 2.** Human MCP-1 percent recovery from different sample types.

<u>Sample</u>	<u>Spike Level (pg/ml)</u>		
	<u>85</u>	<u>268</u>	<u>500</u>
Serum (n=5)	82%	77%	84%
EDTA Plasma (n=5)	79%	85%	88%
Heparin Plasma (n = 5)	108%	100%	101%

**Linearity:** To determine assay linearity, five samples were spiked with a known value of recombinant MCP-1 in serum, EDTA plasma, and heparin plasma. Each sample was serially diluted with standard diluent to produce values that are within the assay range (Table 3).

**Table 3.** Assay linearity from different sample types.

<u>Dilution</u>		<u>Serum</u>	<u>EDTA plasma</u>	<u>Heparin plasma</u>
1:2	Average % of Expected	88	97	91
	Range (%)	82 - 95	93-103	79-99
1:4	Average % of Expected	93	110	115
	Range (%)	84 - 97	104-114	112-120
1:8	Average % of Expected	86	110	90
	Range (%)	83 - 116	106-118	83-97

## Reference

1. *Immunoassay: A Practical Guide*, Chan and Perlstein, Eds., 1987, Academic Press:New York, p71.

This product ("Product") is warranted to operate or perform substantially in conformance with published Product specifications in effect at the time of sale, as set forth in the Product documentation, specifications and/or accompanying package inserts ("Documentation") and to be free from defects in material and workmanship. Unless otherwise expressly authorized in writing, Products are supplied for research use only. No claim of suitability for use in applications regulated by FDA is made. The warranty provided herein is valid only when used by properly trained individuals. Unless otherwise stated in the Documentation, this warranty is limited to one year from date of shipment when the Product is subjected to normal, proper and intended usage. This warranty does not extend to anyone other than the original purchaser of the Product ("Buyer").

**No other warranties, express or implied, are granted, including without limitation, implied warranties of merchantability, fitness for any particular purpose, or non infringement. Buyer's exclusive remedy for non-conforming Products during the warranty period is limited to replacement of or refund for the non-conforming Product(s).**

There is no obligation to replace Products as the result of (i) accident, disaster or event of force majeure, (ii) misuse, fault or negligence of or by Buyer, (iii) use of the Products in a manner for which they were not designed, or (iv) improper storage and handling of the Products.

Current versions of product instructions are available at [www.thermo.com/pierce](http://www.thermo.com/pierce). For a faxed copy, call 800-874-3723 or contact your local distributor.

© 2009 Thermo Fisher Scientific Inc. All rights reserved. Unless otherwise indicated, all trademarks are property of Thermo Fisher Scientific Inc. and its subsidiaries. Printed in the USA.

---

**Data Templates**

	1	2	3	4	5	6	7	8	9	10	11	12
A												
B												
C												
D												
E												
F												
G												
H												

	1	2	3	4	5	6	7	8	9	10	11	12
A												
B												
C												
D												
E												
F												
G												
H												